Temperature Controlled Shaking

Leading Safety Standards

Superior Ease of Use

Reduced Cost of Ownership

Operating Manual

Page 2

Test Tube Shaker Hei-MIX Multi Reax

Operating manual must be read before initial start-up.

Please follow the safety instructions provided.

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About this document	4
Version	4
About this manual	4
Terms and signs used.....	5
Safety Instructions	6
General safety instructions.....	6
EU Declaration of Conformity.....	6
Intended use.....	6
Installation / Electrical safety	7
Personnel qualification.....	7
Operating company's obligations.....	7
Installation site	7
Changes to the unit	8
Safety for the personnel	8
Safety during use	8
Disposal	8
Device Description	9
Device overview	9
Overall view.....	9
Control panel	9
Start-Up	10
Setting up device	10
Set up.....	10
Power cord	10
Connect/disconnect power cord	10
Switch unit on and off	11
Carousel attachment.....	12
Attach vessels	12
Insert vessels.....	13
Operation	15
Operation and controls	15
Continuous shaking.....	16
Timer controlled shaking.....	16
Repetition	16
Edit time during shaking operation	16

Cleaning and Maintenance	17
Cleaning.....	17
Maintenance	17
Troubleshooting	17
Possible errors and how to resolve	17
Assembly	18
Electrical connections.....	18
MULTI-REAX.....	18
Exchange carousel attachment	19
Disassembly and Storage	20
Dismantling, transportation and storage	20
Dismantling	20
Transportation and storage	20
Accessories and Spare Parts	21
Scope of delivery.....	21
Accessories.....	21
Attachments	22
Technical data	22
Service	23
Contact / Technical Service	23
Warranty	24
Confirmation of condition of unit.....	25

Version

Version

This manual describes the function, operation and maintenance of the Hei-MIX Multi Reax test tube shaker.

Version	Alteration date
0.2	03/2017

About this manual

This operating manual is a component of the device described here.

- Please read this manual carefully and obey all safety and warning instructions.
- Ensure that every operator reads this manual.
- Ensure that this manual is accessible for every operator.
- Pass on the operating manual to the subsequent owner.

For the current version of this operating manual in pdf format, please go to:

<http://www.heidolph-instruments.com/Support/Operation Manuals/Shakers & Mixers/Reax and Multi Reax/Multi Reax>

In addition, please observe the regional regulations.

Terms and signs used

In this manual you will find standardized terms and signs to warn you of possible dangers or give important hints. Please take special notice to these signs to avoid personal injury or damage to the unit.

The following terms and type of signs (basic drawing elements) are used:

Symbol	Additional term / Description
Warning signs 	The yellow triangle indicates hazardous situations. It is used in combination with the following terms: DANGER: Indicates a hazardous situation which, if not avoided, will result in irreversible injury or death. WARNING: Indicates a hazardous situation which, if not avoided, could result in serious injury or material damage. CAUTION: Indicates a hazardous situation which, if not avoided, may result in a minor injury or material damage.
Prohibitory signs 	Do not do: The red circle indicates a situation that should be avoided under all circumstances to avoid injury or damage.
Mandatory signs 	Must do: The blue circle indicates important information. Please obey to avoid property damage.

Other signs used

Symbol	Description
→	Handling instruction, action required
✓	Result of action
▪ ▪ ▪	List of information
a. b. c.	List of variants

➤ General safety instructions

EU Declaration of Conformity

This device complies with the following EC-Directives*:

- 2006/42/EC Machinery Directive
- 2014/30/EU Electromagnetic Compatibility Directive

*Also see attachment „EU Declaration of Conformity“.

The device has been constructed according to state-of-the-art technology and recognized safety regulations. However, risks may still arise during installation, operation and maintenance.

→ Please ensure the operating manual is available at all times.

The device may only be used under the following circumstances:

- Only operate the device, if it is in full working order.
- Ensure all operators of the device possess the necessary safety and risk awareness.
- Operate the device according to the instructions stipulated in this manual only.
- If there is something you do not understand, or certain information is missing, ask your manager or contact the manufacturer.
- Do not do anything on the device without authorization.
- Only use the device according to its intended use.

Intended use

The device is intended for use by trained and authorized personnel only.

The device is suitable for the following use:

- mixing
- shaking
- blending of solvents

The operation of the device is suitable in one of the following locations:

In any research laboratory

Improper use

Any use which deviates from the device's intended use is considered to be improper. The manufacturer does not accept liability for any damages resulting from non-permitted uses. The risk is carried by the operator alone.

Installation / Electrical safety

- The device may only be connected when the mains voltage corresponds to the information on the type plate of the unit.
- The mains connection must be accessible at all times.
- Repairs may only be performed by a qualified electrician.
- Never operate the unit with a damaged power cord.
- Always turn the unit OFF and disconnect mains power before performing any maintenance or service.

Personnel qualification

- The device may only be operated by qualified persons.
- The device may only be operated by individuals who have been instructed in its proper use by qualified persons.
- The device may only be operated and maintained by persons who are of legal age.
- Other personnel may only work with the unit under continuous supervision of an experienced qualified operator.
- This manual must be read and understood by all persons working with the device.
- The personnel must receive special safety instructions in order to guarantee responsible and safe work procedures.

➤ Operating company's obligations

Installation site

- The device must be positioned in a suitable location.
- The device must be installed sufficiently stable.
- All screw connections must be securely tightened.
- Do not operate the device in proximity of highly flammable and explosive substances.
- Installation and operation of the device is only permitted in facilities which are fitted with the appropriate laboratory equipment (e.g. with air extraction units).
- We recommend operation in fume hood enclosures.
- The device may only be operated in enclosed spaces and under the following environmental conditions:

Ambient temperature	5 - 31 °C at 80 % relative humidity
	32 - 40 °C decrease linearly till max. 50 % relative humidity
Installation altitude	0 - 2,000 m (6,500 feet) height above sea level
Degree of pollution	2
Overvoltage category	II

Changes to the unit

- No unauthorized changes may be made to the unit.
- No parts may be used which have not been approved by the manufacturer.
- Unauthorized changes result in the EC Declaration of Conformity losing its validity, and the unit may no longer be operated.
- The manufacturer is not liable for any damage, danger or injuries that result from unauthorized changes or from operating the unit other than as described in this manual.

Safety for the personnel

- Ensure that only qualified personnel operate the device.
- Observe the following regulations:
 - Laboratory guidelines
 - Accident prevention regulation
 - Ordinance on Hazardous Substances
 - Other generally accepted rules of safety engineering and occupational health
 - Local regulations

➤ Safety during use

- Wear the appropriate protective clothing when working on the device (clothing, protective glasses and if necessary, safety gloves).
- Do not use the device in potentially explosive areas. The device is not protected against explosion. There is no explosion or ATEX protection available.
- Do not operate or assemble devices in the vicinity which are emission or radiation sources (electromagnetic waves) for the frequency range (3×10¹¹ Hz to 3×10¹⁵ Hz).
- Avoid putting pressure on the display when you are not operating the device.
- Eliminate errors immediately.
- Do not use abrasive material to clean the glass surface. Only wipe with damp cloth.
- Always switch the device off after use.

➤ Disposal

- Check the device components for hazardous substances and solvents.
- Clean all components before disposal.
- Dispose of the device according to the appropriate national legal regulations.
- Dispose of the packaging material in accordance with the appropriate national legal regulations.


➤ Device overview

Overall view

Control panel

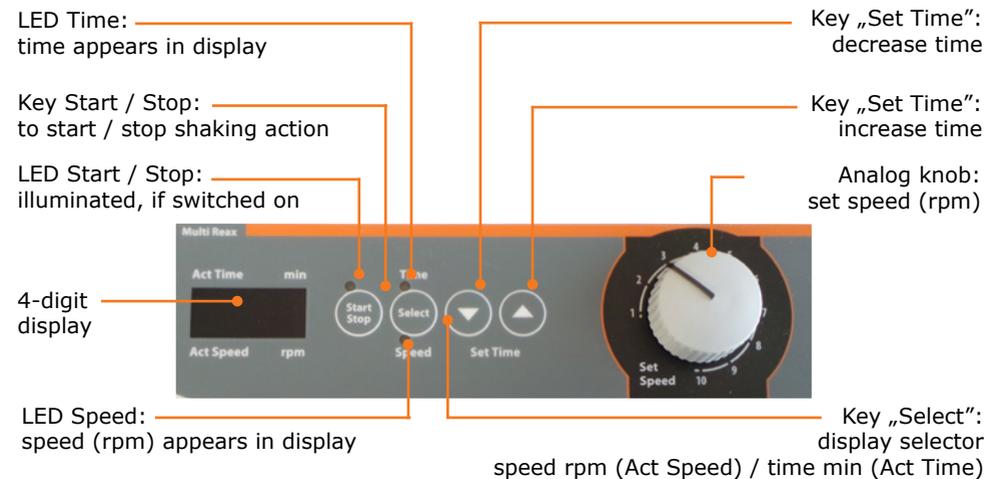➤ Setting up device

Warning: **Risk of slipping device**

As a result of vibrations generated during operation the device might slide off the table top.

The surface must be smooth and clean and should be properly leveled.

Set up

- Locate the shaker on a stable, horizontal surface.
- Clean surface and feet with a damp cloth and ethanol regularly.

➤ Power cord

Connect/disconnect power cord

Connect plug to device

The power cord comes with a three-wire plug and a recess on the bottom side.


Unit plug

Power cord

The socket is located on the back side of the unit.

- Connect plug to the socket with the recess facing down.
- Push plug up against socket until securely attached.
- Connect plug to power socket.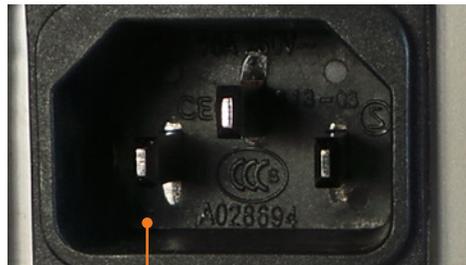

Unit socket

Disconnect power cord

- Disconnect plug from power socket first.
- Then disconnect plug from socket on unit.

➤ Switch unit on and off

The ON/OFF switch is situated on the right side of the device beneath the control panel.

Switch unit on

- Press the ON/OFF switch on the upper side.
- ✓ The unit is switched on.

Switch unit off

- Press the ON/OFF switch on the lower side.
- ✓ The unit powers off.

Switch OFF

Switch ON

➤ Carousel attachment

Attach vessels

Caution:
Risk of breakage and spilling

If the rubber tension ring is damaged your vessels may fall out of the attachment and spill.

Control the rubber tension ring regularly. Replace it eventually.

Vessel sizes

- Carousel attachment for 26 test tubes
 - round vessels, Ø 10 – 16 mm
 - max. length 160 mm
- Carousel attachment for 12 test tubes
 - round vessels, Ø 16 – 32 mm
 - max. length 120 mm.

Attach vessels

- Plug vessels or tubes into holding pockets against preload of a rubber tension ring.
- ✓ This rubber tension ring is going to safely secure the tubes in the holding position.

Carousel attachment

Rubber tension ring

Holding pockets

- To ease installation, lift rubber tension ring by hand before introducing vessels.

Insert vessels



Caution:
Risk of breakage and spillage


If vessels are not inserted properly they may fall out of the attachment and spill.

Refer to recommendations below when inserting vessels.


- Ensure that rubber tension ring holds vessel in center.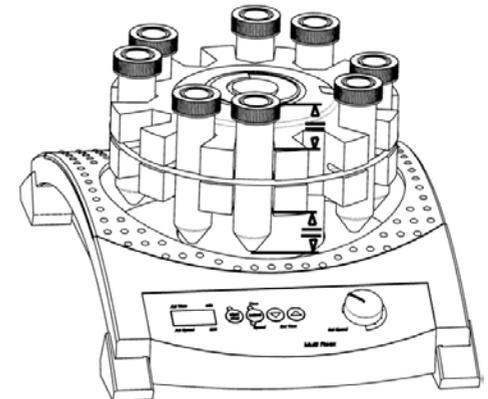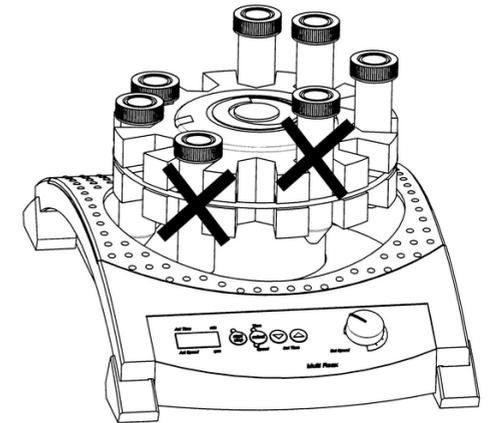

→ Never attach different vessel sizes randomly. The smaller diameter vessel will not be securely fastened!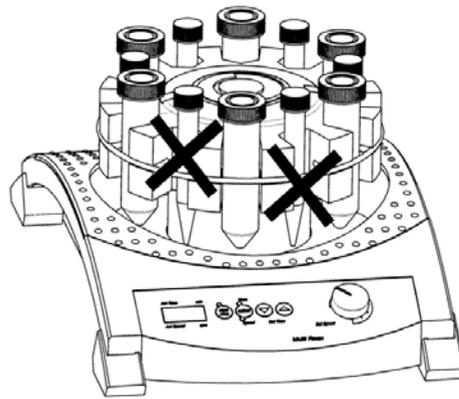

→ Form groups of vessels with same diameter and leave one space between groups open.

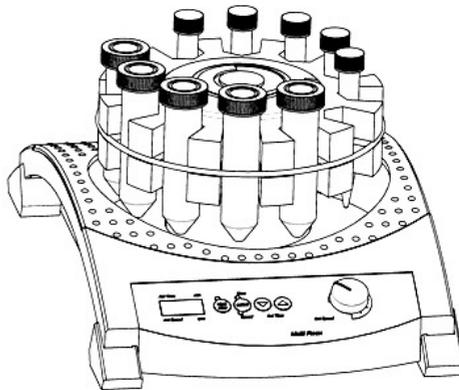➤ Operation and controls

Warning:
Risk of injury!

Returning power after a power cut off will restart shaking mode. There is a risk of injury in the vicinity of moving parts of device. Before working with the device or in the vicinity ensure all controls are switched to "0" or "OFF" mode, when device is not moving.

Caution:
Risk of breakage and spillage

If the vessels are not attached securely they may fall out and spill the sample.

Prior to switching the unit on ensure that all vessels are securely fastened.

Warning:
Risk of poisoning

Open vessels and too high vibration frequencies may result in samples splashing.

Close all vessels tight especially if sample is dangerous or toxic.

Adjust vibration frequency step by step until you have reached the required rpm settings..

Wear safety glasses and especially with dangerous and toxic samples appropriate protective clothing.

Before switching the unit on ensure that no vessel can strike against the housing.

Depending on the quantity and the load of vessels, certain speed levels between 1,200 to 1,500 rpm may create a resonance. Increase or decrease speed by a few rpm until the resonance disappears.

Continuous shaking

- Turn device ON with ON/OFF switch.
- Set time with ↑ and ↓ keys to 000 (depress ↑ and ↓ together).
- Start continuous operation by hitting the Start / Stop key.
- Start and Speed LEDs turn on, shaking frequency is displayed (4-digit display).
- Set shaking frequency with knob.
- Hitting Start / Stop key will stop shaking.

Timer controlled shaking

- Set time between 0 and 999 minutes with Set Time ↑ or Set Time ↓ keys. Short hits on the keys change time setting slowly, whereas depressing keys is going to change time rapidly. Depressing the Set Time ↑ and Set Time ↓ keys resets the timer to 000.
- Depressing the Start / Stop key starts shaking at the frequency selected with knob and for the time interval set before. Start LED turns ON.
- Time lapse appears in a 4-digit display. Time LED turns ON.
- Current shaking action may be discontinued with the Start / Stop key. Actual time won't change. Hitting this Start / Stop key continues shaking at the moment it had been discontinued.
- In case of a power supply failure or interruption (item turned OFF), this factor is lost.
- Hitting the Select key during timer-controlled operation, you may change display from time to shaking frequency and vice versa.
- ✓ Speed LED turns ON. Timer LED flashes to indicate timer operation mode.
- ✓ Timer LED also flashes, when timer-controlled operation was discontinued with the Start / Stop key.

Repetition

Repeating last timer controlled operation does not require new time setting, as long as you did not touch the master switch.

Edit time during shaking operation

- Time setting may be changed during shaking, while timer is running and even when timer controlled operation was discontinued.
- ✓ Memory timer will keep initial value in its memory.

A power cut will stop the unit. All settings will be saved. With power return the device is switched on automatically. Shaking must be started manually.

Cleaning

All surfaces of the unit allow for cleaning with a damp cloth and if required a mild soap lotion.

Attention: Damage to the surfaces

All surfaces can be damaged by improper cleaning.

Never use:

- Chlorine bleach or any cleaning agent including chlorine
- Ammonia
- Abrasive cleaning agents such as cleaning rags, scrubbing agents or any other agents which include metal components

As a result of vibrations generated during operation the device might slide off the table top.

To ensure a stable position clean surface and feet with a damp cloth and ethanol regularly.

Maintenance

The unit is maintenance-free. Any necessary repair must be performed by an authorized Heidolph distributor.

Please contact Heidolph Instruments or your local Heidolph Instruments distributor.

Possible errors and how to resolve

Malfunctions	Possible reason	Troubleshooting
Unit won't shake	No power	<ul style="list-style-type: none"> ▪ Check size and shape of plug and compatibility to your electrical socket. ▪ Check circuit breakers
	Light diode defective	Contact your local Heidolph Instruments distributor
Unit stops shaking	Pre-set timer elapsed	Restart timer, if necessary
	Thermal motor circuit breaker triggered by motor overload.	Wait 20 minutes, decrease overall load
	<ul style="list-style-type: none"> ▪ Mechanical parts broken ▪ Electronic failure 	Contact your local Heidolph Instruments distributor

If you experience a malfunction which can not be resolved, please contact your authorized Heidolph Instruments distributor immediately.

Electrical connections

- Repairs of any kind are allowed by **qualified professional electricians only**. Any improper repair can result in a dangerous situation. Contact your local Heidolph Instruments distributor for any repair you may have.

Warning:
Electric shock!

This unit must only be connected to a grounded electrical socket.

The unit and voltage must match. The type label on the back side of the unit provides all voltage specifics.

When shipped the unit is grounded. In case the original plug is changed, the new plug must have a protective conductor!

MULTI-REAX

Rating plate

Color code for electrical connections:

Europe		North America	
GREEN/YELLOW	PE: Protective conductor (Earth)	GREEN	PE: Protective conductor (Earth)
BLUE	N: Neutral conductor	WHITE	N: Neutral conductor
BROWN	P: Phase	BLACK	P: Phase

The unit is connected with the power cord supplied to the electrical socket. It plugs into the housing on the rear side of the unit.

For countries utilizing a plug other than the standard one supplied:

- The plug supplied may only be changed by a professional electrician.
- If utilizing an adaptor ensure that it is grounded and approved by the local regulations.

Exchange carousel attachment

The item arrives with the carousel attachment for 26 test tubes installed. You may replace it with the carousel attachment for 12 test tubes at ease.


- Remove all vessels from carousel attachment.
- Remove carousel attachment by pulling with both hands.
- Plug new carousel attachment in place.

Check for location of the two "noses". Carefully slip carousel attachment over these "noses". Be sure not to damage the part. These "noses" secure the carousel on the shaker.

Check rubber tension ring for proper seat. This ring secures your vessels in place. Seat ring in slot of carousel on condition.

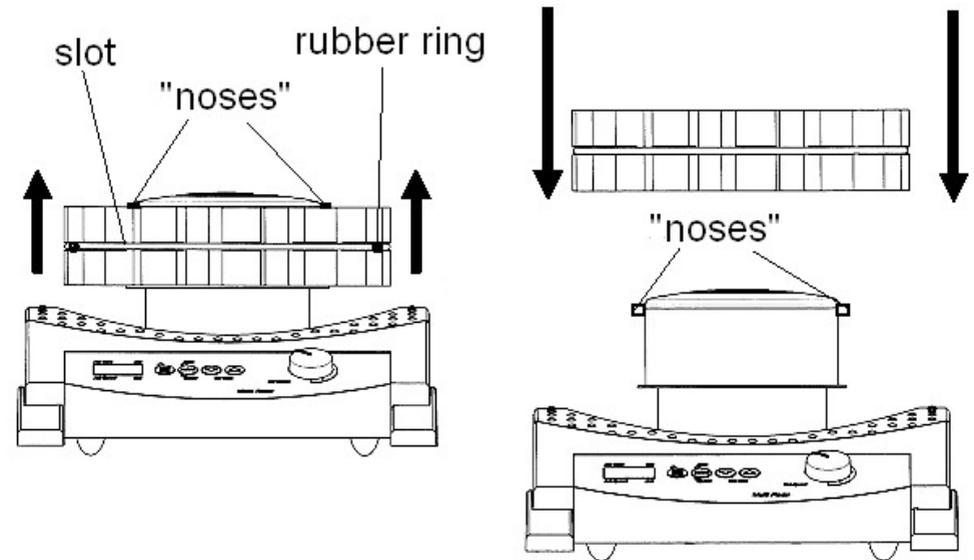➤ Dismantling, transportation and storage

Dismantling

**Warning:
Risk of injury!**

While the unit is plugged in you might be at harm if the device is accidentally switched to ON mode.

Prior to dismantling switch the unit off and pull the plug from the electric socket. Ensure that the unit is not switched on again.

- ✓ The unit is switched off.
- Disconnect plug from power socket.
- Disconnect plug from socket on the unit.

Transportation and storage

- Store and transport the unit and its components only if they were emptied and cleaned,
- Store and transport the unit and parts of unit in the original packing material or alternatively in an adequate container to prevent damages.
- Seal the packing carefully against unauthorized or mistaken opening.
- Store the unit in a dry and frost-free place.

Improper storage and transportation may result in damages to the system and the mechanical components.

Avoid any kind of shocks during transportation.

➤ Scope of delivery

Component	Quantity	Product number
Hei-MIX Multi Reax	1	545-10000-00
Power cord	1	14-007-003-81
Carousel attachment for 26 test tubes, 10 - 16mm Ø	1	549-01600-00
Carousel attachment for 12 test tubes, 16 - 32mm Ø	1	549-03000-00
Operating Manual English/German	1	01-005-004-28
Warranty registration / Confirmation of condition	1	01-006-002-78
EC Declaration of Conformity	1	01-001-025-06

➤ Accessories

Component	Quantity	Product number
Carousel attachment for 12 test tubes, 16 - 32 mm Ø	1	549-03000-00
Rubber tension ring for carousel attachment 26 tubes	1	23-08-06-01-75
Rubber tension ring for carousel attachment 12 tubes	1	23-08-06-01-76

Additional accessories are shown in the general catalog or on our website at

- www.heidolph-instruments.com

➤ Technical data

Multi Reax

Standard supply voltage	230 V/50 Hz or 115 V/60 Hz
Input power (W)	50
Protection class (DIN EN 60529)	IP 30
Drive	asynchronous motor
Overheat protection	self-resetting
Rotation speed (rpm)	150 - 2,000
Rotation speed setting	digital
Motion	circular vibrating
Orbit (mm)	3
Load capacity (kg)	1.5
Operating mode	automatic or continuous
Weight (kg)	9.8
Dimensions (W x h x d)	270 x 410 x 172
Permissible ambient temperature (C°)	<ul style="list-style-type: none"> ▪ 5 - 40 °C ▪ 80 % relative humidity, no condensation at 5 - 31 °C, linear decreasing to 50 % at 40 °C

➤ Contact / Technical Service

Questions / Repair work

If any aspect of installation, operation or maintenance remains unanswered in the present manual, please get in touch with the following address.

For repairs please call Heidolph Instruments or your local authorized Heidolph Instruments distributor.

Warning: Danger of poisoning!

Contaminated units can lead to severe injury or death of our employees!

When shipping items for repair that may have been contaminated by hazardous substances, please

- advise exact substance
- take adequate protective action towards our parts receiving and service personnel
- mark the pack in accordance with Ordinance on Hazardous Substances

Included at the end of this manual you will find a „Confirmation of condition of unit“.

→ Kindly copy and fill in this form and submit it prior to shipping the unit for repair.

Contact details:

Heidolph Instruments Germany

Heidolph Instruments GmbH & Co. KG
Technical Service
Walpersdorfer Str. 12
D-91126 Schwabach / Germany

Tel.: +49 - 9122 - 9920-74
Fax: +49 - 9122 - 9920-84

E-Mail: service@heidolph.de
www.heidolph.de

Heidolph Instruments North America

Phone: 1-866-650-9604
E-mail: service@heidolph.com
www.heidolphNA.com

Heidolph Instruments United Kingdom

Phone: 01799 - 5133-20
E-mail: service@radleys.co.uk
www.heidolph-instruments.co.uk

All other countries

You will find contact details of your local Heidolph distributor at www.heidolph.com

➤ Warranty

Heidolph Instruments provides a 3-year warranty for the products described here (excluding glass and wearing parts) if you register using the warranty card enclosed or online (www.heidolph.com). The warranty is valid from the point of registration. The serial number is also valid without registering.

The warranty covers part and manufacturing defects.

In the case of a part or manufacturing defect, the device shall be repaired or replaced free of charge under the terms of the warranty.

Heidolph Instruments shall not assume liability for any damages incurred as a result of improper handling or transport.

Warranty claim?

➔ Please inform Heidolph Instruments or your local Heidolph distributor should you wish to make a warranty claim.

➤ Confirmation of condition of unit

➔ In the case of repair, copy and complete the Confirmation of condition of unit and send it to Heidolph Instruments.

1. Details about the unit

Product number _____
 Serial number _____
 Reason for repair _____

2. Has the device been cleaned, decontaminated/sterilized?

Yes _____ No _____

3. Is the unit in a condition which does not represent any health threats for the staff of our service department?

Yes _____ No _____

If not, which substances has the unit come into contact with?

4. Legally binding declaration

The customer is aware of being legally liable to Heidolph Instruments for any damages arising from incomplete and incorrect information.

Date _____ Signature _____

Company stamp

Please note

The shipper is responsible for the return of the goods in well packed condition, suitable for the mode of transport.

Sender information

Name, first name _____
 Company _____
 Department, research group _____
 Street _____
 Zip code, city _____
 Country _____
 Phone _____
 E-mail _____

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This is a translation of the original operating manual. We reserve the right to make any technical modifications without prior notification. This documentation is not subject to revision; the latest version of this documentation can be found on our homepage. The copyright for all texts and images is held by Heidolph Instruments GmbH & Co. KG.
© Heidolph Instruments GmbH & Co. KG